

ZP400 Capsule & Tablet Counting Machine



Instructions:

ZP400 Capsule & Tablet Counting Machine is used for counting the drugs or foods like the capsule, tablet, grain, etc. It is mainly used in the fields such as pharmaceutical, hospital and foodstuff. The whole machine is made of the stainless steel. The counting plate adopts unique design; it can count and fill the tablet and capsule. It adopts “Eccentric screen” structure and frequency converter, overcoming the problem of loud noise, easy damaged, difficult adjusted, etc. It is characterized by small volume, light weight, accurate counting, simple operation, etc. The contact surface of countered material is also made of stainless steel. It is the ideal counting equipment for vial, bag or can products in various relevant industries.

Technical Parameters:

Model	ZP400
Counting Plate Diameter	400 mm
Hole Quantity Of The Plate	5-100 (Optional)
Production Capacity	12-26 bottles(bags)/ min
Power	220V/50Hz/0.55 Kw
Packing Size	600×500×700 mm
Packing Weight	55 Kg

Manual

1. Then a good power supply, so that the number of tablets of disc counter-clockwise rotation.

SINOPHAM CO., LTD

ADD: NO.128, NORTH ZHANGJIABANG ROAD, CHINA(SHANGHAI) PILOT FREE TRADE ZONE, SHANGHAI, CHINA

TEL: 0086-170-9142-7771

MAIL: MARKETING@SINOPHAM.COM

HTTP: WWW.SINOPHAM.COM

2. Into an appropriate capsule, open the power supply, will be transferred to a suitable voltage switch position, so that electromagnetic vibration to reach the right intensity, so that the number of tablets capsules disk hole easy access to a few tablets, while dynamic capsule with a brush, so that the number of tablets of hole both capsules.

3. Bottling, it will bottle feed into the capsule out of the mouth or the mouth of the bottle with your fingers gently touch the micro-switch, the number of tablets of disc rotation to all fall into the capsule under the feed mouth, and then immediately release micro-switch.

4. Into the next count work, using brush dynamic capsule

Caution:

1. The number of tablets of the height of the disk should be the head capsule length of the same or slightly higher than the number of tablets of Capsule disk, not too low or too high, otherwise turn will ruin capsule is too low can be higher than the pad with the Calibrating the middle part and to the middle hole ended up with the shaft after the match abandoned DVD films, is too high please lathe to a few tablets of the bottom plate washer (do not be confused too low).

2. Collar disk on the number of particles must be adjusted to the appropriate height, and bottom should be the head capsule to maintain a distance of about 1.5-2mm, too low, hampering the capsules will be too high.

3. The number of tablets not all covered with holes in the capsule not to be pressed micro-switch, otherwise the position will go to Circlip broke capsule. Counting all the holes in the capsule is not available when covered with a brush dynamic capsule, so that the number of tablets are capsule hole.

4. Damp after the deformation of the capsule can not count on the machine.

5. The number of tablets not pressure plate, or not working properly; lift the machinery can not be carried in a few tablets of the plate, otherwise it would loose a few tablets of disc damage.

6. Capsule to be locked together completely, the length of a unified, without fail capsules.

Note:

1. to lift the machinery must not carry on in a few tablets of disc hopper, or cause damage to the machine!

2. on the machine feet: You can install the 4 screws and tighten the nut after the feet as the machines used.